

E. coli Water Test Kit Instructions

Use the pipette to procure water sample.

Measure 5 ml of the water sample into a graduated cup.

Pour the water sample into the media.

Reseal the media bottle and shake.

Uncap the tube and remove the hardening stick; place on a clean, dry surface.

Pour the sample/media into the tube.

Insert the hardening stick into the tube.

Cap the tube; be sure the cap is tight.

Insert the tube in incubator box/stand, upright.

Allow the media to solidify (15-20 minutes).

The tube may be placed inside the box for incubation Allow 24 hours for culture.